

Deep UV Extension for the MicroTime 200



- Excitation at 266 nm with compact diode based laser
- Quartz lenses and UV compatible optics
- UV objective and UV detector for optimal performance
- Parallel UV and VIS detection
- Label-free Fluorescence Lifetime Imaging Microscopy (FLIM)
- Label-free protein detection through intrinsic fluorescence

VIS excitation:

375 nm, 405 nm, ..., 640 nm




Deep UV excitation:

- 266 nm
- < 100 ps pulse width
- variable repetition rates from 1 MHz to 80 MHz



Modified main optical unit
with quartz optics, for parallel
VIS and UV detection

to microscope

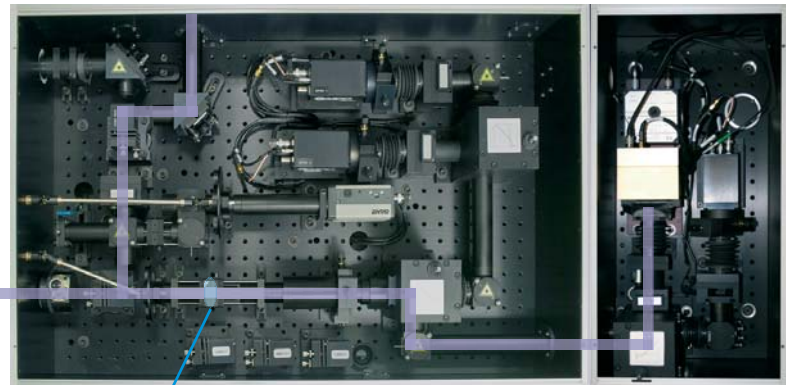

Modified microscope body:
quartz objective
(glycerine immersion)




Quartz tube lens

Detection extension unit with optimized
PMT for UV detection:
42 % detection
efficiency at 350 nm


Deep UV Extension for the MicroTime 200

Applications

FLIM with streptavidin-coated beads
(\varnothing 500 nm, 1 streptavidin = 24 tryptophans)
immobilized on quartz coverslips

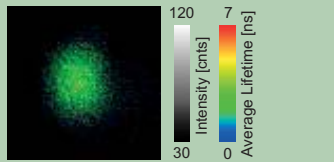


1.5 x 1.5 μ m

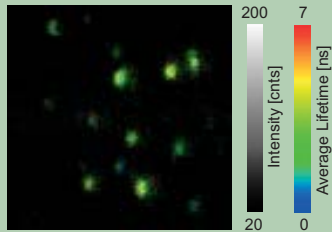


13 x 13 μ m

**Label-free FLIM
in the deep UV**

Label-free FLIM with biological cells
→ aromatic amino acids within the proteins
(mainly tryptophans) become visible

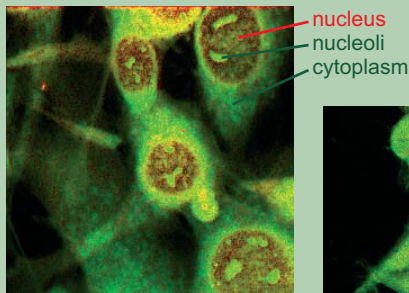


3T3-L1 cells, fixed (methanol)
80 x 80 μ m

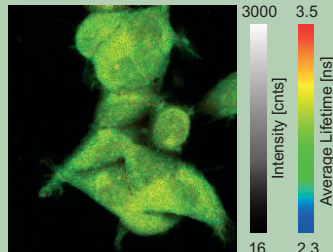

HEK293 cells, fixed
(paraformaldehyde)
73 x 79 μ m

*Sample courtesy of Astrid Tannert,
University of Leipzig, Germany*

**Dyes with high quantum efficiency
in the UV: laser dyes as single
emitters**

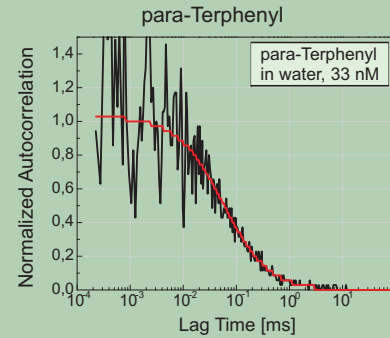


- diffusion time: $t(\text{diff}) = 57 \mu$ s
- molecular brightness: 510 Hz per molecule
- data acquisition time: approx. 10 min

**Benchmark:
FCS in the deep UV**

Streptavidin-coated beads
(\varnothing 120 nm)

approx. 1000 streptavidin per bead
(1 streptavidin = 24 tryptophans)

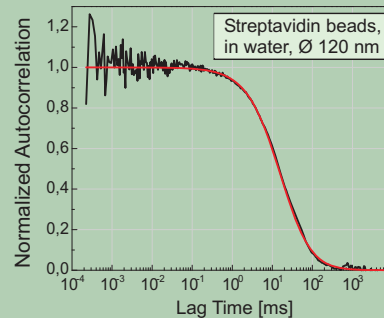

- diffusion time: $t(\text{diff}) = 16$ ms
- molecular brightness: 1.2 kHz per molecule
- data acquisition time: approx. 5 min

- excitation: 266 nm (PicoQuant), 20 MHz, < 100 ps pulse width
- quartz objective, 40x, NA 0.6, glycerine
- optical filters: Z266RDC, 300 nm longpass
- detection: PMT, 42 % detection efficiency at 350 nm

All Information given here is reliable to our best knowledge. However, no responsibility is assumed for possible inaccuracies or omissions. Specifications and external appearances are subject to change without notice. Trademarks or corporate names are used for explanation and identification, to the owner's benefit and without intent to infringe.

© PicoQuant GmbH, June 2014



PicoQuant GmbH
Rudower Chaussee 29 (IGZ)
12489 Berlin
Germany

Phone +49-(0)30-6392-6929
Telefax +49-(0)30-6392-6561
Email info@picoquant.com
WWW <http://www.picoquant.com>